Image Quant™ TL GxP 10.2

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Windows 版

2023.06

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The screenshot shows the Cytiva software interface. On the left, there is a 'FILTERS' sidebar with various status filters. The main 'Projects' area displays a list of projects, with '001 test - ACTIVE' selected and highlighted in blue. Below it are other projects like 'test 5 - COMPLETE', 'test 2 - COMPLETE', 'test 3 - QC-CHECKED-OUT', 'test 1 - COMPLETE-CHECKED-OUT', and 'test - COMPLETE-CHECKED-OUT'. On the right, the '001 test' details are shown, including a 'HISTORY' section with a list of events and their timestamps. At the bottom right, the 'ACTIONS' section contains buttons for 'Check Out', 'Check Out (R/O)', and 'Request Signoff', with the latter being highlighted by a red rectangle. Below the actions is an 'AUDIT LOG' section with a 'Report' button.

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The screenshot shows a 'Notes' dialog box. It has a text input field containing the text 'signoff request'. Below the input field are two buttons: 'OK' and 'Cancel'. The 'OK' button is highlighted with a red rectangle.

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The screenshot shows the Cytiva software interface after the signoff request. The '001 test' project is now highlighted in blue and its status is 'AWAITING-SIGNOFF', indicated by a red exclamation mark icon. A red arrow points to the project name. The 'HISTORY' section on the right now includes a new entry: 'Request Signoff demo2 signoff request' with a timestamp of '17 May 2022 16:00'. The 'ACTIONS' section now includes a 'Cancel Signoff Req.' button. The 'AUDIT LOG' section remains the same.

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
The screenshot displays the Cytiva software interface. On the left, there is a 'FILTERS' sidebar with various filters like 'Last 7 days', 'Active', and 'All'. The main area is titled 'Projects' and lists several projects. The first project, '001 test - AWAITING-SIGNOFF', is highlighted with a blue border and a red arrow pointing to it. Below it are other projects like 'test 5 - COMPLETE', 'test 2 - COMPLETE', 'test 3 - QC-CHECKED-OUT', 'test 1 - COMPLETE-CHECKED-OUT', and 'test - COMPLETE-CHECKED-OUT'. On the right, there is a '001 test' sidebar showing a 'HISTORY' of actions and an 'ACTIONS' section with buttons for 'Check Out (R/O)', 'Cancel Signoff Req.', and 'Sign Off'. The 'Sign Off' button is highlighted with a red box.

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
This is a close-up view of the 'ACTIONS' section from the previous screenshot. It shows three buttons: 'Check Out (R/O)', 'Cancel Signoff Req.', and 'Sign Off'. The 'Sign Off' button is highlighted with a red rectangular box. Below the actions section is an 'AUDIT LOG' section with a 'Report' button.

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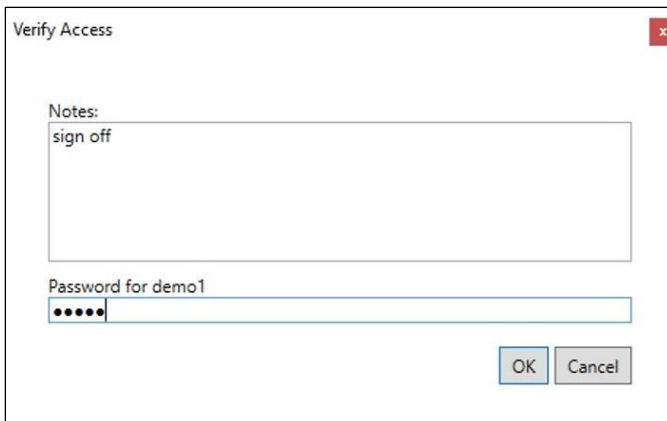The image shows a 'Verify Access' dialog box. It has a title bar with a close button (X). Inside, there is a 'Notes:' label followed by a text input field containing the text 'sign off'. Below this is a 'Password for demo1' label followed by a password input field with six dots. At the bottom right, there are two buttons: 'OK' and 'Cancel'.

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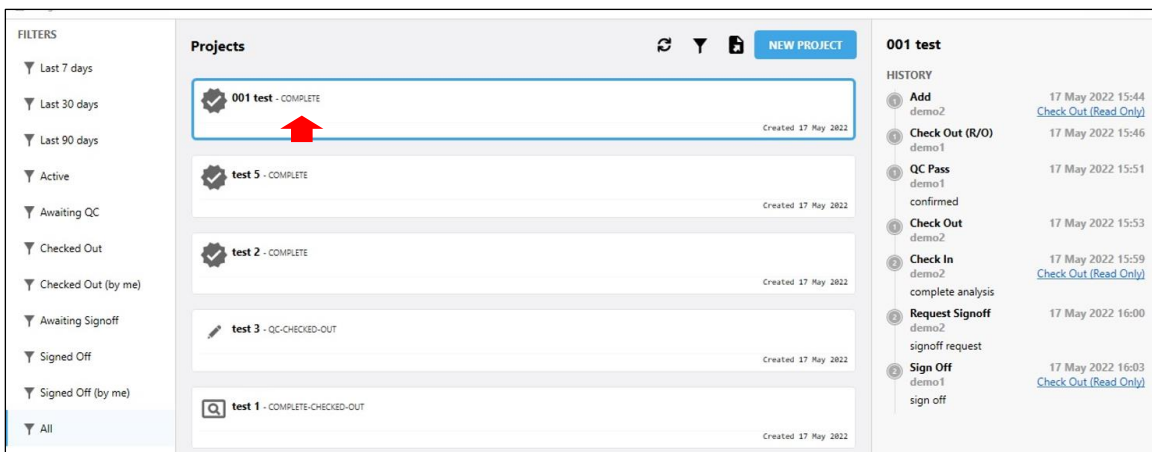
The image is a screenshot of a web application interface. On the left is a 'FILTERS' sidebar with various status filters like 'Last 7 days', 'Active', etc. The main area is titled 'Projects' and shows a list of project cards. The top card is '001 test - COMPLETE' and is highlighted with a blue border and a red arrow pointing to it. To the right of the project list is a 'HISTORY' section for the selected project, showing a list of actions with timestamps and links like 'Check Out (Read Only)'. A 'NEW PROJECT' button is visible in the top right of the main area.

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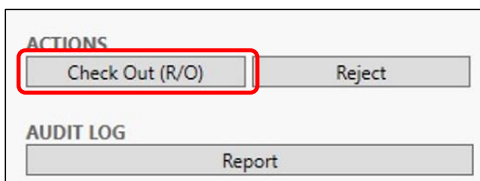
The image shows a close-up of the 'ACTIONS' menu. There are two buttons: 'Check Out (R/O)' and 'Reject'. The 'Check Out (R/O)' button is highlighted with a red rectangular box. Below the actions menu is an 'AUDIT LOG' section with a 'Report' button.

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
The image is a screenshot of the 'Projects' page, similar to the one above. The project '001 test - COMPLETE-CHECKED-OUT' is now highlighted with a blue border and a red arrow pointing to it. The 'AUDIT LOG' section is not visible in this view.

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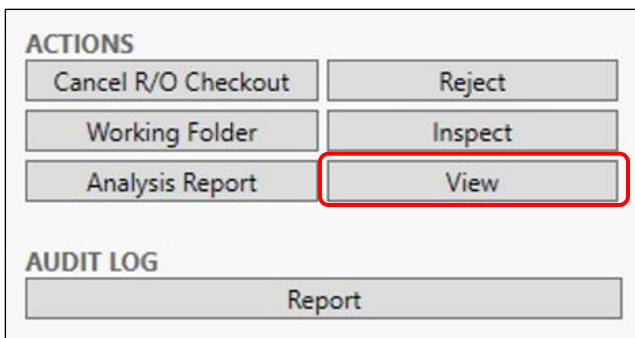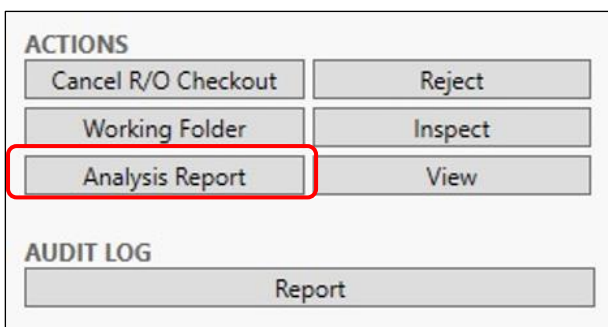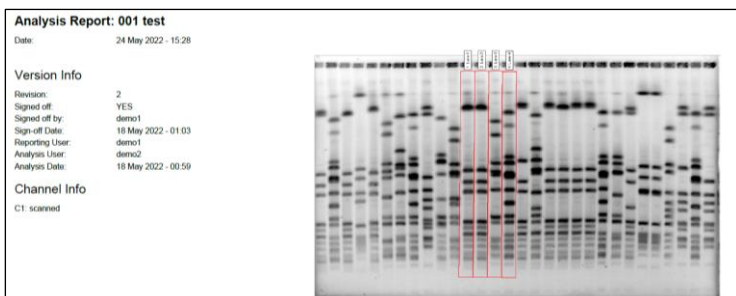
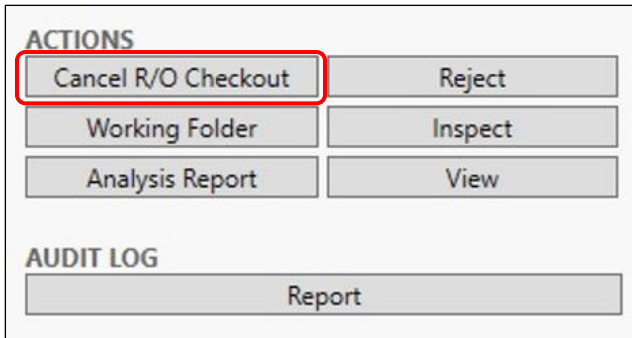
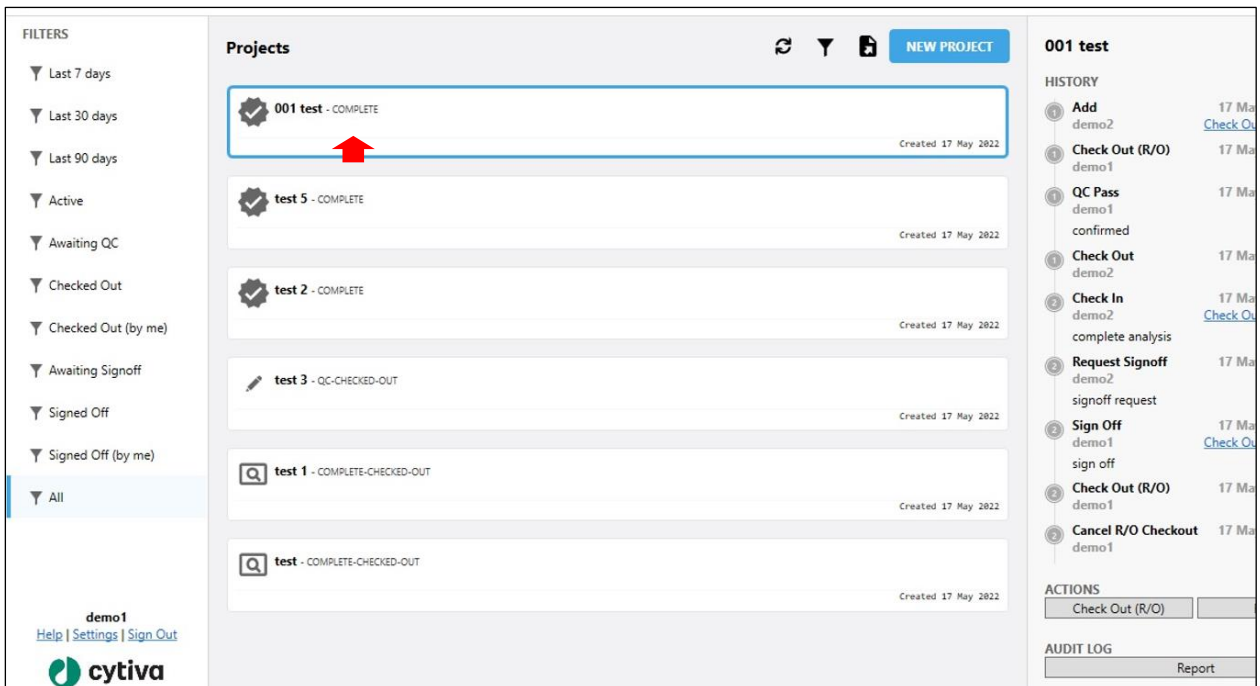
The screenshot shows the Cytiva software interface. On the left, there is a 'FILTERS' sidebar with various status filters. The main area displays a list of 'Projects' with details like '001 test - COMPLETE' and 'Created 17 May 2022'. On the right, the '001 test' details are shown, including a 'HISTORY' section with a list of events and their timestamps. At the bottom right, there is an 'AUDIT LOG' section with a 'Report' button highlighted by a red box.

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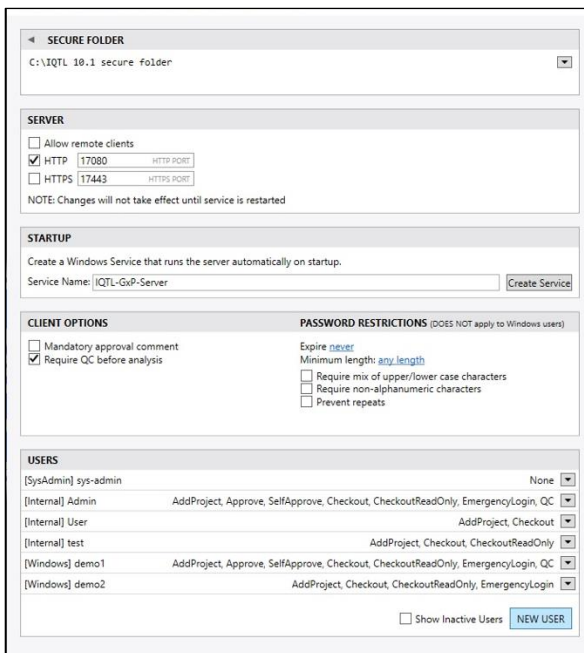
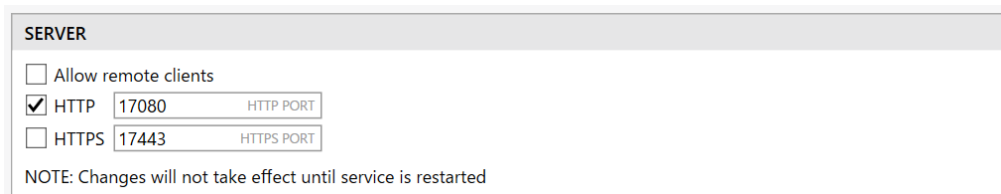
CLIENT OPTIONS	PASSWORD RESTRICTIONS (DOES NOT apply to Windows users)
<input type="checkbox"/> Mandatory approval comment <input type="checkbox"/> Require QC before analysis	Expire never Minimum length: any length <input type="checkbox"/> Require mix of upper/lower case characters <input type="checkbox"/> Require non-alphanumeric characters <input type="checkbox"/> Prevent repeats

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5. USERS

USERS	
[SysAdmin] sys-admin	None ▼
[Internal] Admin	AddProject, Approve, SelfApprove, Checkout, CheckoutReadOnly, EmergencyLogin, QC ▼
[Internal] User	AddProject, Checkout ▼
[Internal] test	AddProject, Checkout, CheckoutReadOnly ▼
[Windows] demo1	AddProject, Approve, SelfApprove, Checkout, CheckoutReadOnly, EmergencyLogin, QC ▼
[Windows] demo2	AddProject, Checkout, CheckoutReadOnly, EmergencyLogin ▼
<input type="checkbox"/> Show Inactive Users NEW USER	

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7 IQTL GxP regulatory requirements

Regulatory authorities require pharmaceutical development and manufacturing companies to use quality-assured equipment and to follow detailed and carefully documented analytical and manufacturing procedures. The purpose of this control effort is to ensure consistent and reliable quality in pharmaceutical products that reach the market. GxP regulations describes requirements for companies and laboratories, and include guidelines on how standard protocols can be used for safe electronic record keeping. IQTL GxP has functions that supports compliance to the **21 CFR Part 11** regulation by Food & Drug Administration (FDA) and **EU GMP Annex 11**.

The following functions support IQTL GxP 21 CFR Part 11:

- **Audit trails (21 CFR 11.10 (e))**
 - Project audit history – complete log of user activity
 - Analysis log – detailed application history
 - Access log – full list of all successful / failed login attempts
- **Limiting system access (21 CFR 11.10 (d))**
 - Robust password access
 - Ability to use Windows User Access Control for authentication
 - Fine grained feature access based on permissions
 - Client/server approach prevents user access to secure folder
- **Electronic signatures (21 CFR 11, Subpart B)**
 - Approval functionality limited to configured users
- **Password restrictions (21 CFR 11, 11.200 (a))**
 - Full compliance with passwords complexity requirements
- **Emergency logins (21 CFR 11, 11.200 (3))**
 - Two users (with appropriate permissions) can provide credentials to pose as another for emergency situations
- **Version History for change control (21 CFR 11, 11.10 (k))**
 - Resilient data integrity enforced throughout the lifetime of a project

- Detailed log of all operations performed on a project

Note: IQTL GxP is intended to address GxP requirements in compliance with 21 CFR part 11 and EU GMP Annex 11. It is the user organization's responsibility to make sure that the software is appropriately validated within the context of the user's particular environment, and to make sure that the other Part 11 or EU GMP Annex 11 compliance elements are implemented appropriately.

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate.

ImageQuant TL is a trademark of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

Windows, Active Directory, and Excel are registered trademarks of Microsoft Corporation. Mac and OS X are trademarks of Apple Incorporated. Acrobat is a trademark of Adobe Systems Incorporated.

All other third-party trademarks are the property of their respective owners.

© 2021 Cytiva

Any use of ImageQuant TL is subject to Cytiva Standard Software End-User License Agreement for Life Sciences Software Products. A copy of this Standard Software End-User License Agreement is available on request.

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.

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