Image Quant TL

(ver.10.0)

Gel and Blot Analysis 簡易操作手順

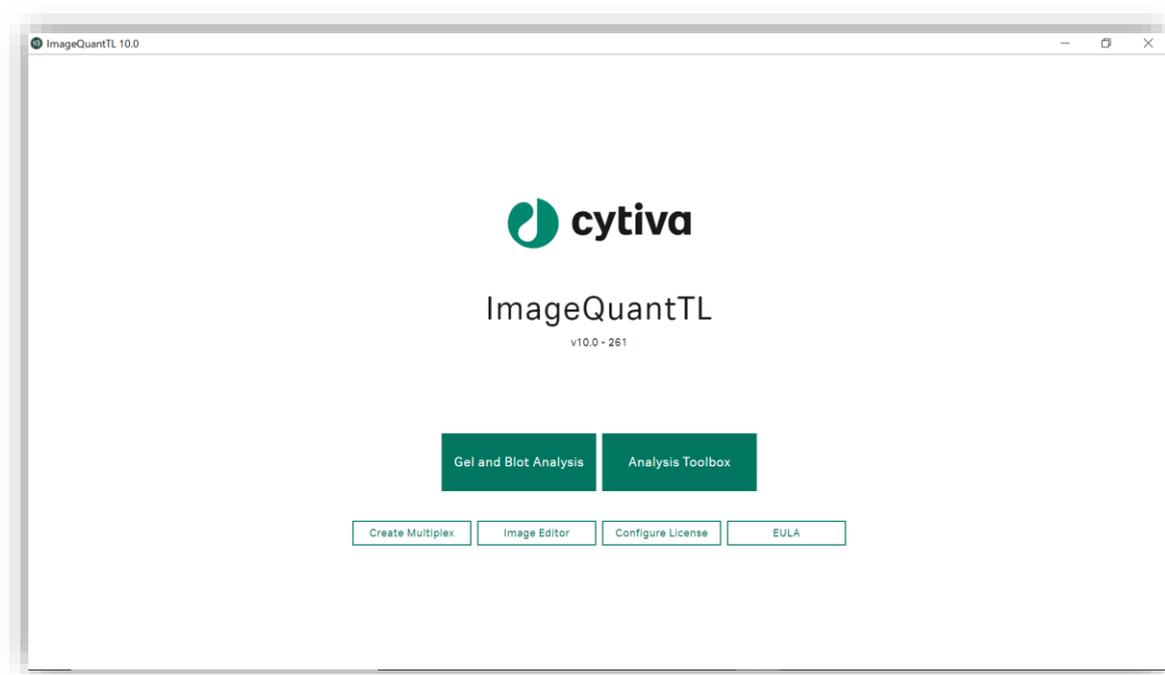
Image Quant TL 操作手順書

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Normalization Factor = $\frac{\text{Volume of largest overall HK band in reference channel}}{\text{Volume of HK band in each lane in reference channel}}$ > or = 1.0

Normalized Volume = Normalization Factor x Volume of each lane in active channel

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Normalization Factor = $\frac{\text{Volume of most abundant lane in reference channel}}{\text{Volume of each lane in reference channel}}$ > or = 1.0


Normalized Volume = Normalization Factor x Volume of each lane in active channel

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Zoom Image Fit Display [Multiple Channels]

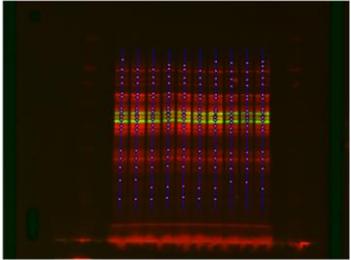Image Lanes Bands MW Compare Normalize

Lane	1	2	3	4	5	6	7	8	9	10
Normalized Band Volume	0.000000	0.000000	0.000000	0.000000	0.000000	0.000000	0.000000	0.000000	0.000000	0.000000

Define Normalization

Normalize band areas across your gel

Method: Total Protein

Reference Channel: 20180831 3 color, manual probe

Normalization Factor Denominator: Total Band Volume

For each lane, a normalization factor NF is calculated by dividing the total volume of the most abundant lane by its own total volume. Measured band volumes in the reference channel. Total volume is either the volume of the whole lane, or the sum of all band volumes for the lane. Bands are assigned normalized values by multiplying their measured volume by the NF for the lane.

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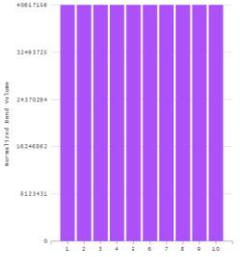
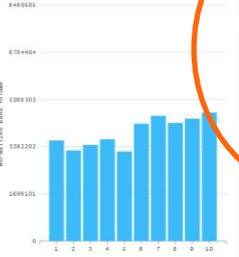
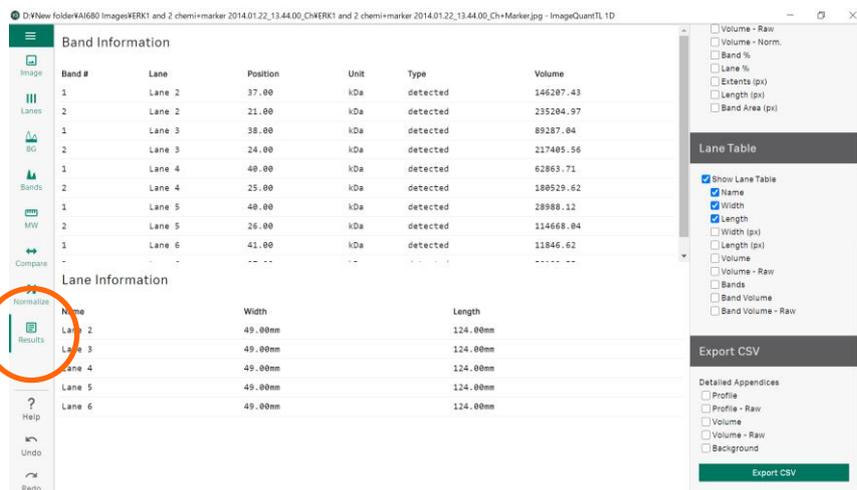
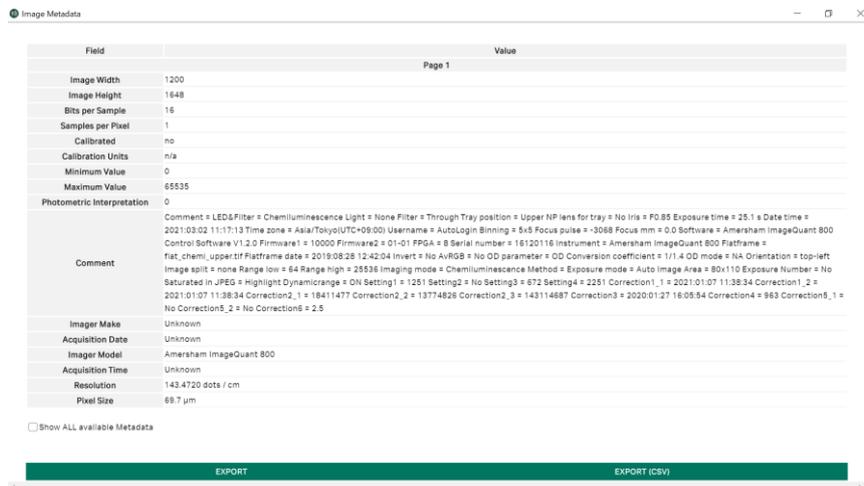
Field	Value
Page 1	
Image Width	1200
Image Height	1648
Bits per Sample	16
Samples per Pixel	1
Calibrated	no
Calibration Units	n/a
Minimum Value	0
Maximum Value	65535
Photometric Interpretation	0
Comment	Comment = LED&Filter = Chemiluminescence Light = None Filter = Through Tray position = Upper NP lens for tray = No Iris = F0.85 Exposure time = 25.1 s Date time = 2021:03:02 11:17:13 Time zone = Asia/Tokyo(UTC+09:00) Username = AutoLogin Binning = 5x5 Focus pulse = -3088 Focus mm = 0.0 Software = Amersham ImageQuant 800 Control Software V1.2.0 Firmware1 = 10000 Firmware2 = 01-01 FPGA = 8 Serial number = 16120116 Instrument = Amersham ImageQuant 800 Flatframe = Flat_chemi_upper.tif Flatframe date = 2019:08:28 12:42:04 Invert = No Av/RGB = No OD parameter = OD Conversion coefficient = 1/1.4 OD mode = NA Orientation = top-left Image split = none Range low = 64 Range high = 25536 Imaging mode = Chemiluminescence Method = Exposure mode = Auto Image Area = 80x110 Exposure Number = No Saturated in JPEG = Highlight Dynamicrange = ON Setting1 = 1251 Setting2 = No Setting3 = 672 Setting4 = 2251 Correction1_1 = 2021:01:07 11:38:34 Correction1_2 = 2021:01:07 11:38:34 Correction2_1 = 18411477 Correction2_2 = 13774826 Correction2_3 = 143114887 Correction3 = 2020:01:27 16:05:54 Correction4 = 963 Correction5_1 = No Correction5_2 = No Correction6 = 2.5
Imager Make	Unknown
Acquisition Date	Unknown
Imager Model	Amersham ImageQuant 800
Acquisition Time	Unknown
Resolution	143.4720 dots / cm
Pixel Size	69.7 µm

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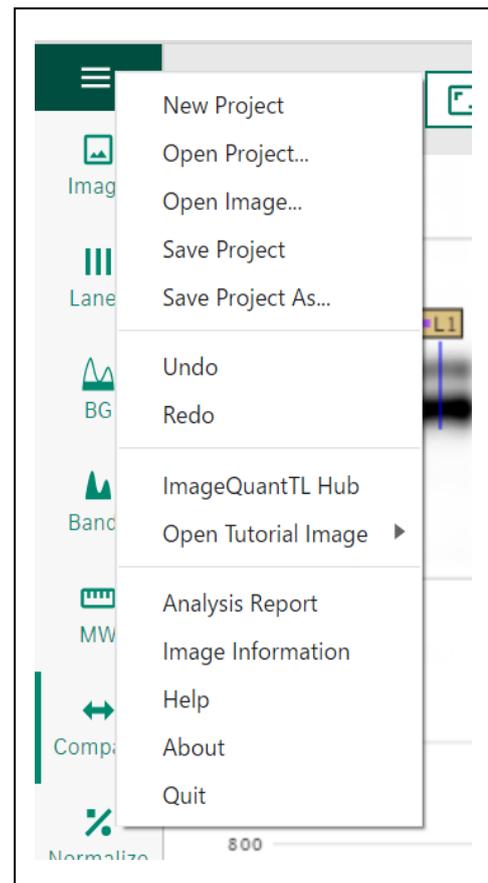
Trademarks

For local office contact information, visit cytiva.com/contact

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate.

Amersham and ImageQuant are trademarks of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

Apple is a trademark is a trademark of Apple Incorporated. Windows is a trademark of Microsoft Corporation.

All other third-party trademarks are the property of their respective owners.

© 2021 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request.

Contact your local Cytiva representative for the most current information.

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